


Gel Extraction Cutting Tool

With the Gel Extraction Cutting Tool, DNA and RNA bands can be accurately excised from agarose gels. For one-handed, safe and easy gel cutting, this tool is the disposable alternative to other cutting tools such as razor blades and scalpels without the risk of cross contamination.




IMPORTANT BEFORE USE!


1. **DO NOT** squeeze the tool before dispensing the gel in step 3.
2. This tool was designed for 1 time use only. Discard the tool after use.

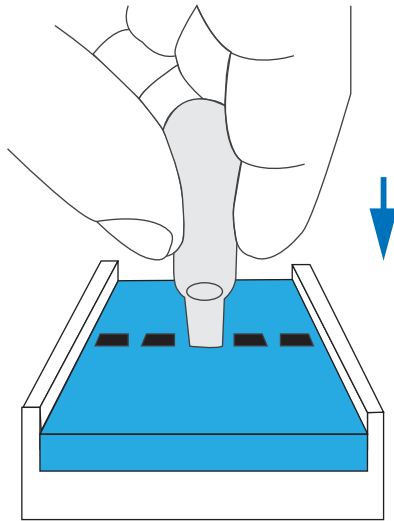

1. Plunge cutter firmly into gel
DO NOT SQUEEZE

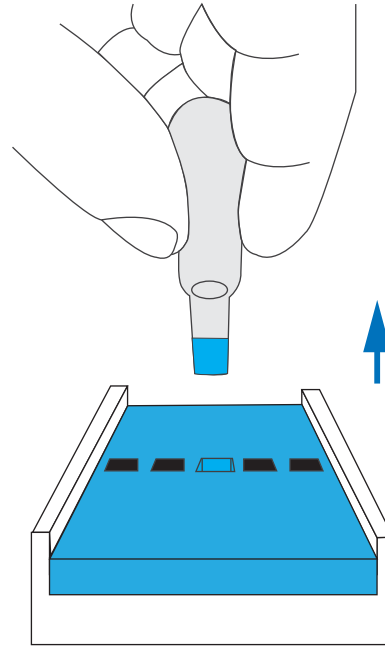

2. Remove selected gel slice
DO NOT SQUEEZE

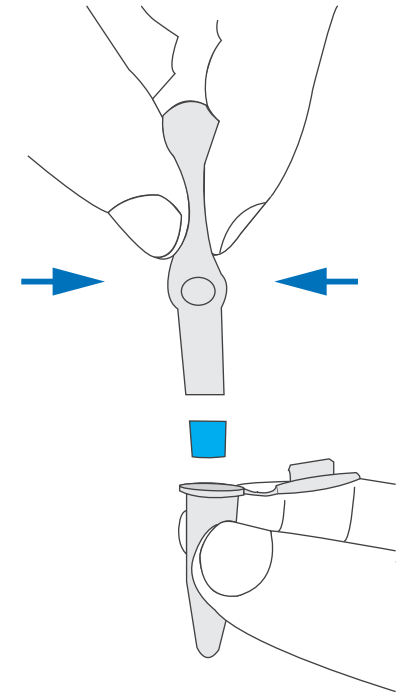

3. Squeeze quickly and firmly to dispense gel slice